

See more with Optical Z-sectioning

A module for the NEW IN Cell Analyzer 1000

The Optical Z-sectioning Module is a combination of hardware and proprietary software control that removes out-of-focus light above and below the focal plane. Suitable for use in both end-point and kinetic versions of the new IN Cell Analyzer

1000, the Optical Z-sectioning Module allows the generation of high-quality 2D and 3D images of fluorescently labeled cellular and tissue samples for analysis where high background is problematic.



Fig 1. IN Cell Analyzer 1000 is a modular and fast analysis platform for cell-based assays.



Enhanced sample flexibility

Optical Z-sectioning allows publication-quality imaging due to improved signal-to-noise ratio and increased contrast for samples such as fluorescent tissue samples, cell aggregates (e.g. stem cells), and some migration assays. Optical Z-sectioning also eliminates the need for wash steps in some cellular assays: decreasing time to result and increasing productivity.

Subcellular localization and 3D reconstruction

Exportable image formats from an IN Cell Analyzer 1000 enabled with the Optical Z-sectioning Module allow high-quality volume reconstructions to be made. Specific cellular events can also be clearly localized.

User-selectable modes for increased choice

The new IN Cell Analyzer 1000 enabled with optical Z-sectioning can be switched easily between widefield and Z-sectioning acquisition modes during operation to allow the most productive use of the instrument.

Flexible choice of fluors


IN Cell Analyzer 1000's lamp-based illumination maximizes the choice of fluors that can be imaged in widefield or Z-sectioning mode. Up to four wavelengths can be acquired sequentially enabling a wide range of assays and multiplexed experiments to be performed.

Build the instrument you need


The Optical Z-sectioning Module is an optional upgrade for both endpoint and kinetic versions of the new IN Cell Analyzer 1000. For maximum versatility in live- and fixed-cell assay development, other upgradeable modules are available including Temperature Control, Liquid Handling, Microscope Slide Imaging, and Transmitted Light Imaging.

A wide range of cell imaging applications

Automated multiwavelength image acquisition and analysis can be performed in both widefield and optical Z-sectioning modes in a wide range of assay types including tissue imaging, tissue assays, migration assays, and organelle and protein trafficking studies.

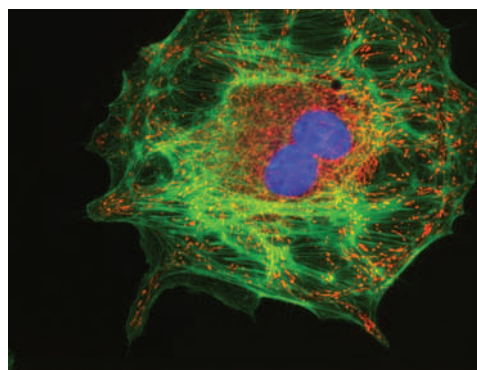


A. Widefield

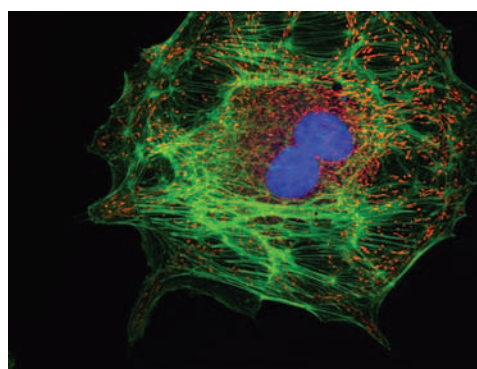


B. Optical Z-section

Fig 2. FluoCells™ prepared slide contains bovine pulmonary artery endothelial cells (BPAEC). The mitochondria are stained with MitoTracker™ Red CMXRos, F-actin is labeled with BODIPY™ FL phalloidin, and the nuclei are labeled with DAPI.

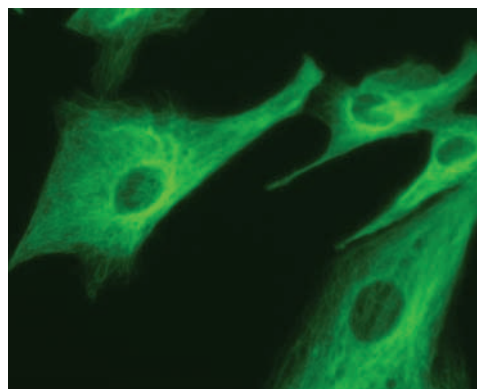

A. Widefield

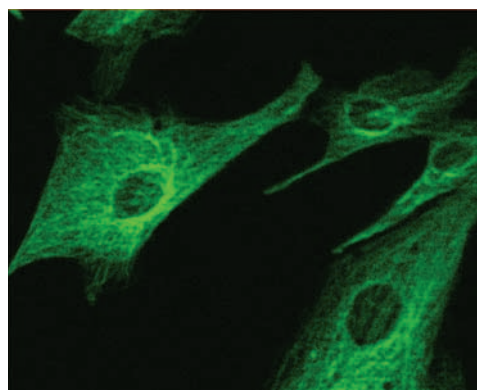

B. Optical Z-section

Fig 3. Bovine pulmonary artery endothelial cells stained with mouse monoclonal anti-tubulin antibody to label microtubules (FITC).

General system specifications		
Camera	Pixel size	6.45 × 6.45 μm
	Field of view at 10×	0.603 mm ²
	Cooling	-30 °C
	Pixel layout	1392 × 1040
Focus	Method	Confocal laser sensor
	Focus time	< 400 ms
	Objective positioner resolution (Piezo objective positioner)	0.2 μm
Illumination	Epifluorescent lamp	100 W xenon
	Transmitted light	LED
Magnifications	4×/0.20 plan apo	
	10×/0.45 plan apo	
	20×/0.45 ELWD plan fluor	
	40×/0.60 ELWD plan fluor	
Plate positioning (XY stage)	Resolution	+/- 5 mm
	Speed (well-to-well speed on a 96-well microplate)	< 800 ms
Filter selection	Excitation filters	6
	Emission filters	6
Plate compatibility	SBS standard footprint microplates including 6-, 12-, 24-, 48-, 96-, and 384-well plates. Glass slides	
Computer	Standard Windows XP Professional (SP2) PC	
Kinetic system specifications		
Reagent dispensing	Configuration	Single needle dispenser
	Plate compatibility	Specified 6-, 12-, 24-, 48-, 96-, and 384-well plates
	Addition/transfer volume	10–100 μl
Temperature control	Minimum setting	Ambient + 5 °C
	Maximum setting	42 °C
Miscellaneous specifications		
Electrical	Voltage	110–240 V
	Frequency	50–60 Hz
Instrument physical specifications	Depth	970 mm
	Width	670 mm
	Height	710 mm
	Weight	~70 kg
Service cabinet	Depth	610 mm
	Width	560 mm
	Height	660 mm
	Weight	~ 50 kg
Automation	Robotics ready and compatible with all major commercially available laboratory robotic systems	

Ordering information

Product	Code number
Optical Z-sectioning Module for the NEW IN Cell Analyzer 1000*	28-4051-30
IN Cell Analyzer 1000 end-point instrument	28-4051-28
IN Cell Analyzer 1000 kinetic instrument†	28-4051-29
Temperature Control Module	28-4063-95
Liquid Handling Module	28-4066-59
Microscope Slide Imaging Module	25-8098-19
Transmitted Light Module	28-4077-07

Product	Code number
IN Cell Investigator Software, single seat license (1 machine license)‡	25-4089-71
IN Cell Investigator Software, 5 seat licenses (5 concurrent licenses)‡	25-4089-72
IN Cell Dual Stand Monitor	28-4051-31

* The Optical Z-sectioning Module is only compatible with the new IN Cell Analyzer 1000 end-point and kinetic instruments.

† Includes Liquid Handling and Temperature Control Modules.

‡ The IN Cell Investigator Software suite provides a comprehensive solution to high-content image and data analysis by combining the latest version of IN Cell Developer Toolbox and IN Cell Analysis Modules with Spotfire™ DecisionSite™ visualization software. More information can be found in the data file *IN Cell Investigator high-content analysis software v1.0*, GE Healthcare, 28-9009-46, Edition AA (2006).

Asia Pacific	Tel: +85 65 62751830	Fax: +85 65 62751829
Australasia	Tel: +61 2 8820 8299	Fax: +61 2 8820 8200
Austria	Tel: 01 /57606 1613	Fax: 01 /57606 1614
Belgium	Tel: 0800 73 890	Fax: 02 416 8206
Canada	Tel: 1 800 463 5800	Fax: 1 800 567 1008
Central & East Europe	Tel: +43 1 972 720	Fax: +43 1 972 722 750
Denmark	Tel: +45 70 25 24 50	Fax: +45 45 16 2424
Eire	Tel: 1 800 709992	Fax: +44 1494 542010
Finland & Baltics	Tel: +358 9 512 3940	Fax: +358 9 512 39439
France	Tel: 01 69 35 67 00	Fax: 01 69 41 98 77
Germany	Tel: 0800 9080 711	Fax: 0800 9080 712
Greater China	Tel: +852 2100 6300	Fax: +852 2100 6338
Italy	Tel: 02 26001 320	Fax: 02 26001 399

Japan	Tel: 81 3 5331 9336	Fax: 81 3 5331 9370
Korea	Tel: 82 2 6201 3700	Fax: 82 2 6201 3803
Latin America	Tel: +55 11 3933 7300	Fax: +55 11 3933 7304
Middle East & Africa	Tel: +30 210 96 00 687	Fax: +30 210 96 00 693
Netherlands	Tel: 0800-82 82 82 1	Fax: 0800-82 82 82 4
Norway	Tel: +47 815 65 777	Fax: +47 815 65 666
Portugal	Tel: 21 417 7035	Fax: 21 417 3184
Russia, CIS & NIS	Tel: +7 495 956 5177	Fax: +7 495 956 5176
Spain	Tel: 902 11 72 65	Fax: 935 94 49 65
Sweden	Tel: 018 612 1900	Fax: 018 612 1910
Switzerland	Tel: 0848 8028 10	Fax: 0848 8028 11
UK	Tel: 0800 515 313	Fax: 0800 616 927
USA	Tel: +1 800 526 3593	Fax: +1 877 295 8102

www.gehealthcare.com

GE Healthcare Limited
Amersham Place
Little Chalfont
Buckinghamshire
HP7 9NA
UK

GE Healthcare reserves the right, subject to any regulatory approval if required, to make changes in specifications and features shown herein, or discontinue the product described at any time without notice or obligation. Contact your GE Healthcare representative for the most current information. © 2006 General Electric Company - All rights reserved. GE and GE Monogram are trademarks of General Electric Company. FluoCells, Bodipy, and MitoTracker are trademarks of Molecular Probes Inc. The IN Cell Analyzer 1000 and associated analysis modules are sold under license from Cellomics Inc. US patent Nos 6573039, 5989835, 6671624, 6416959, 6727071, 6716588, 6620591 6759206; Canadian patent No 2328194, 2362117, 2,282,658; Australian patent No 730100 European patent 1155304 and other pending and foreign patent applications.

