

Assembly Instructions

1. Insert chiller tubes through the side of the containment chamber.
2. Place the base of the ring stand and stir plate inside the chamber.
 - a. Make sure to have the power cord run through the slot cut in the non-hinged side of the chamber.
 - b. Have the metal poll and electrode plate attached to the base of the stir plate.
3. Feed the electrode wires through the top of the chamber and insert into the electrode plate. **Make sure to properly connect the electrodes to the power supply (Platinum connected to the positive)**
4. Attach the alligator clips and electrodes to the electrode wires.
5. Place the chiller bath onto the stir plate and attach the chillers tubes to the barbed pipe fittings. **Make sure to have the outlet tube from the chiller on the BOTTOM pipe fitting. The INLET tube of the chiller on the TOP pipe fitting.**